

MRC-Holland Technical Support


Portal > Knowledgebase > Coffalyser.Net > Working with Coffalyser.Net > Comparative Analysis > How can I export all sample reports at once?


How can I export all sample reports at once?

This article was retrieved from support.mlpa.com on Friday, 15th February 2019.

This article is part of the [Coffalyser.Net Tips & Tricks](#) collection.

Coffalyser.Net conveniently allows you to export sample reports of all samples simultaneously.

- In the comparative analysis tab, right-click on a sample and select Open Sample Results.
- You can now start the export of all sample reports in three ways:
 1. Navigate to the Sample result overview or Sample report tab, right-click and select Create PDF Reports All Samples.
 2. In any tab, press the keyboard shortcut Ctrl + Alt + R.
 3. In any tab, click the  button in the toolbar.
- In the dialog box that opens, enter a name for the folder in which the reports will be saved in the File name field, and select the location where you want this folder to be created.
- Click Save to create the folder and sample reports.

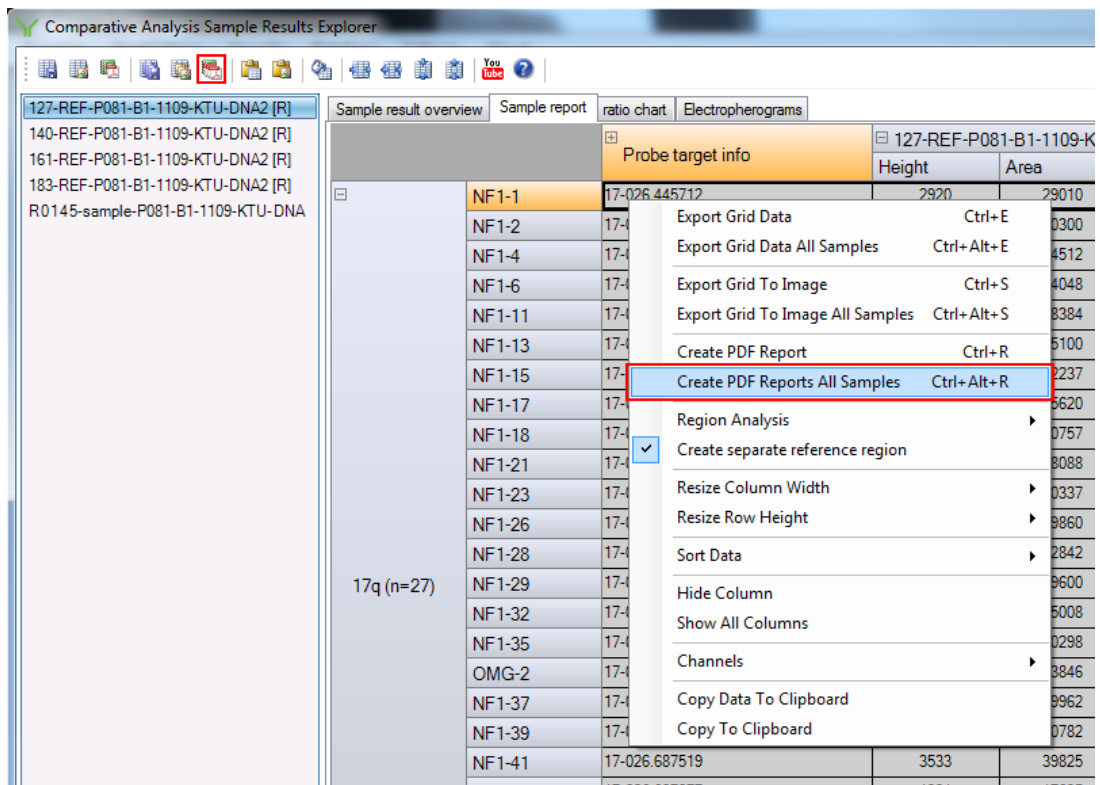

The screenshot shows the 'Comparative Analysis Sample Results Explorer' window. The 'Sample report' tab is active, displaying a table of sample results. A right-click context menu is open over the table, with the option 'Create PDF Reports All Samples Ctrl+Alt+R' highlighted in red. The table contains the following data:

Sample ID	Probe target info	Height	Area
NF1-1	17-026.445712	2920	29010
NF1-2	17-026.687519	3533	39825
NF1-4	17-026.687519	3533	39825
NF1-6	17-026.687519	3533	39825
NF1-11	17-026.687519	3533	39825
NF1-13	17-026.687519	3533	39825
NF1-15	17-026.687519	3533	39825
NF1-17	17-026.687519	3533	39825
NF1-18	17-026.687519	3533	39825
NF1-21	17-026.687519	3533	39825
NF1-23	17-026.687519	3533	39825
NF1-26	17-026.687519	3533	39825
NF1-28	17-026.687519	3533	39825
NF1-29	17-026.687519	3533	39825
NF1-32	17-026.687519	3533	39825
NF1-35	17-026.687519	3533	39825
OMG-2	17-026.687519	3533	39825
NF1-37	17-026.687519	3533	39825
NF1-39	17-026.687519	3533	39825
NF1-41	17-026.687519	3533	39825

Tags

Coffalyser.Net

Disclaimer

The information provided in this material is correct for the majority of our MLPA products. However, for certain applications, the instructions for use may differ. In the event of conflicting information, the relevant instructions for use take precedence.