

# MRC-Holland Technical Support

Portal > Knowledgebase > Coffalyser.Net > Working with Coffalyser.Net > Comparative Analysis > How can I open the results of an individual sample directly from the comparative analysis experiment explorer?

## How can I open the results of an individual sample directly from the comparative analysis experiment explorer?

This article was retrieved from support.mlpa.com on Saturday, 16th February 2019.

This article is part of the [Coffalyser.Net Tips & Tricks](#) collection.

Coffalyser.Net allows you to open the results of individual samples directly from the comparative analysis experiment explorer. It is not necessary to close the experiment explorer first.

To open a sample's results, right-click on the column of a specific sample in the grid and select Open Sample Results. Please note that you cannot right-click on the sample name itself.

		Probe target...	AllSamples	ReferenceSamples	140-I
FRSS (n=7)		n/a	100%	100%	
CAS (n=5)	FMRS	n/a	100%	95%	
	PSLP - Relati...	n/a	OK	OK	
	FSLP - Relati...	n/a	OK	OK	
	RSQ - Refere...	n/a	OK	OK	
	RPQ - Refere...	n/a	Warning	OK	
NF1-1	17-026.445712				
NF1-2	17-026.507125				
NF1-4	17-026.514364				
NF1-6	17-026.532562				
NF1-11	17-026.552886				
NF1-13	17-026.565642				

Tags  
Coffalyser.Net

### Disclaimer

The information provided in this material is correct for the majority of our MLPA products. However, for certain applications, the instructions for use may differ. In the event of

*conflicting information, the relevant instructions for use take precedence.*